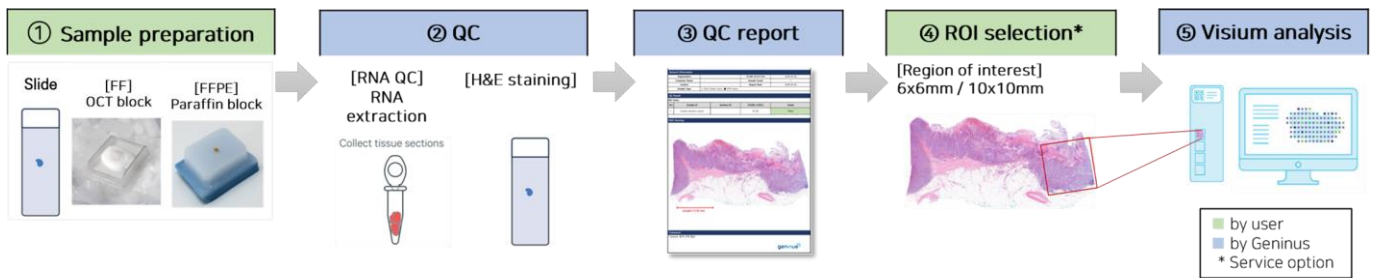
1. Sample preparation protocol for Visium spatial transcriptome analysis (p.2 ~ 5)

2. Sample preparation protocol for Xenium spatial transcriptome analysis (p.6)

For spatial transcriptome analysis			
Visium 6.5 x 6.5 mm	FF OCT block	Embedded in mold & freeze with dry ice	-80 °C (dry ice O)
	FFPE Paraffin block	Embedded in paraffin & store at 4 °C	4 °C (ice pack O)
Visium CytAssist 6.5 x 6.5 mm 10.5 x 10.5 mm	FF Slide (Section: 10 μm)	Slide size: 75 mm x 25 mm x 1 mm A total of 5 slides are required	-80 °C : store the slides in 50 ml conical tube (dry ice O)
	FFPE Slide (Section: 5 μm)	<ul style="list-style-type: none"> ▶ 3 slides for QC (H&E staining + RNA QC + Extra) ▶ 2 slides for main (2 Positively Charged slides: Main + Extra) 	4 °C : store the slides in 50 ml conical tube (ice pack O)
Xenium 10 x 10 mm 10 x 22 mm	FF OCT block	Tissue embedding (OCT) & freeze with dry ice	-80 °C (dry ice O)
	FFPE Paraffin block	Tissue embedding (Paraffin) & store at 4 °C	4 °C (ice pack O)

1. Sample preparation protocol for Visium CytAssist spatial transcriptome analysis

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Table 1. Positively charged slides validated for Visium CytAssist

Item	Length (mm)	Width (mm)	Thickness (mm)
EpreDia Shandon ColorFrost Plus	75.0	25.0	1.0
Fisher SuperFrost/ColorFrost	75.0	25.0	1.0
Sigma-Aldrich Poly Prep Slides	75.0	25.0	1.0
VWR SuperFrost Plus	75.0	25.0	1.0

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*DV200 : Percentage of RNA fragments > 200 nucleotides

*RIN : RNA Integrity Number

1. Sample preparation protocol for Visium CytAssist spatial transcriptome analysis

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a. Incubation using a Section Dryer Oven

- ① Place the slides in a slide drying rack.
- ② Close the lid when incubating the slide in the oven.

b. Incubation using a Thermocycler

- ① Position a Low Profile Thermocycler Adapter on a thermal cycler that is set at the incubation temperature.
- ② Ensure that the Low Profile Thermocycler Adapter is in contact with the thermal cycler surface uniformly.
- ③ When incubating a slide, position the slide on the Low Profile Thermocycler Adapter with the side with tissue facing up.
- ④ Ensure that the entire bottom surface of the slide is in contact with Low Profile Thermocycler Adapter.
- ⑤ DO NOT close the thermal cycler lid when incubating the slide.
- ⑥ Allow Low Profile Thermocycler Adapter to cool down before removing from thermal cycler.

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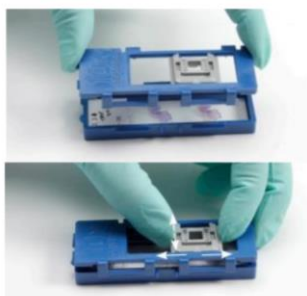
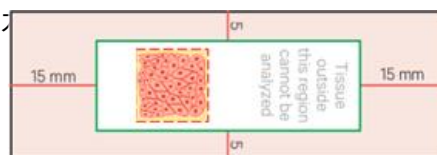
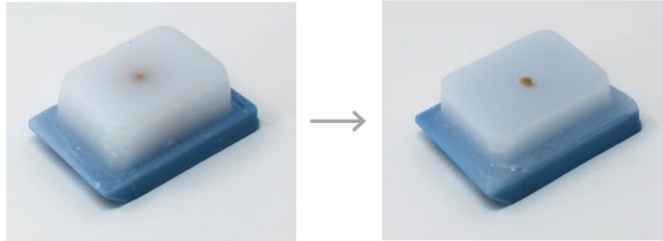
Video : <https://www.youtube.com/watch?v=Wiq96lFyu6s>

3. Sample preparation

3.3 Formalin-fixed paraffin-embedded block


- (1) Collect sample and fix in 10 % formalin or 4 % PFA
- (2) Process the tissue as following (Ethanol → Xylene → Wax)
- (3) Proceed paraffin embedding
- (4) Store at 4 °C or send with ice pack

Exposing the Tissue Block


2. Sample preparation protocol for Xenium spatial transcriptome analysis

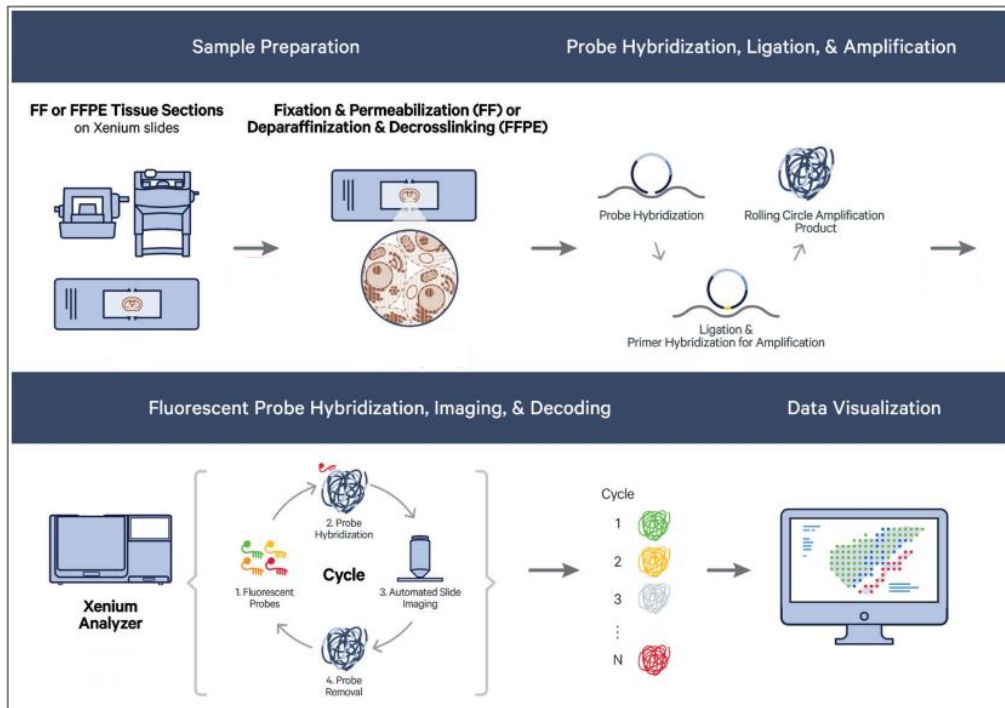1. Service process overview

2. Slide information

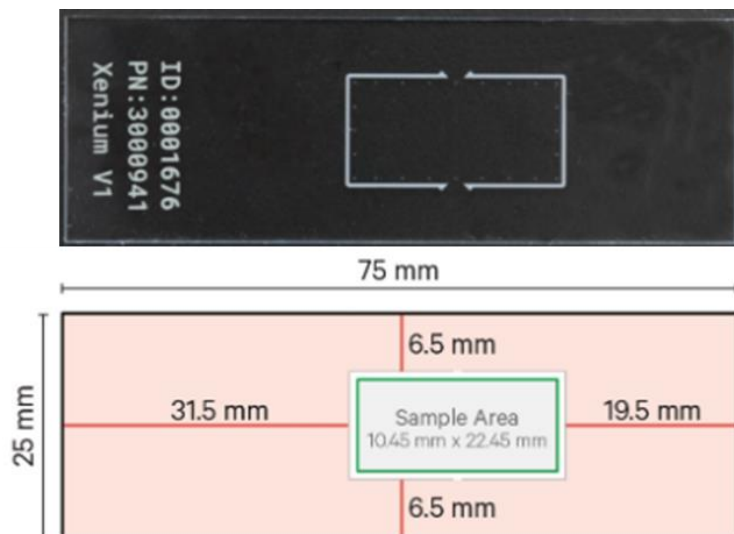- Slide area 10.45 x 22.45 mm
- ROI must be within the green colored area

3. Sample preparation

- FF OCT block & FFPE paraffin block are prepared as above (see Visium CytAssist sample preparation p.4~5)